

## מקלונים למשטחים לשימוש עם צלחות Compact dry

### Compact Dry Swab for surfaces, meat and dry areas

- Unscrew the orange swab and wipe a defined surface area.
- Put the swab back into the tube and screw tightly. Shake the tube and invert to mix.
- Unscrew the cap of the tip. By pressing the middle of the tube squeeze all solution (1ml) on the Compact Dry plates.
- If necessary prepare further series of 10 times sample dilution, and inoculate 1ml each in the middle of Compact Dry plates.

